



您只需一支经济实用的拓普森液体控制器来实现从0.1毫升到100毫升的玻璃和塑料移液管的移液操作

疏水性滤膜防止液体阀门系统渗透,提供提供标准配置的过滤器。拓普森能够121°C的高温消毒。

The economical pipette controller which limits overaspiration. You only need one larger volume pipette Controller for the entire range of bulb and graduated pipettes made of glass and plastic from 0.1 to 100 ml.

Penetration by liquids is effectively reduced by the hydrophobic membrane filter. A replacement filter is provided as standard. The entire macro can be autoclaved at 121°C.



耐用而可靠的大容量Topscien移液器

Durable and reliable large capacity Topscien remove liquid device

Topscien将以往的基本功能理念与先进移液设计和技术融合在一起,使其拥有高品质的外观,同时兼顾移液器的安全性,舒适性和高效性。

The basic function of Topscien will ever ideas and advanced design and technology moving fluid together, make its have high quality appearance, for both the security of moving fluid, comfort and efficiency.



吸液/Filling

挤压橡皮球,可一次性吸入50毫升液体。吸液途径可根据吸八拉到杆的位置连续变化
Squeeze the suction bellows. It allows aspiration up to 50ml with one compression. The fill rate is continuously variable with the pipetting lever.

排液/Delivery

将控制杆向下扳,则液体流出,阀门系统的特殊设计使液体流动并不影响移液精变。
To discharge the liquid, simply press down the lever. The valve system is designed so that flow rate does not affect accuracy.

吹出/Blow-out function

当使用吹出型液管一旦半液型液面停顿不能流出时,按下吹出管内液体。
When using blow-out pipettes, press the convenient blow-out bulb once the meniscus has come to rest in the jet. Pipetting has never been easier.

性能和特点/Performance and feature

- 无需保养设计/Unnecessary maintenance
- 优良的耐化学腐蚀性/Chemical resistance
- 符合人体工程操作舒适/Human engineering design
- 自重轻,结构坚固/Lightweight and Structure strong
- 操作轻便省力/Lightweight and dexterous for operation
- 适合宽范围的移液分配/A wide range volume for liquid controlling

大容量移液器/Large volume pipettes :

产品货号/Cat.No.	产品编号/Code No.	量程/Volume	颜色/Colour
801107	TL01B	0.1ml-100ml	蓝色部件+灰色球型手握球 Blue parts + Gray ball
801108	TL01Y	0.1ml-100ml	黄色部件+灰色球型手握球 Yellow parts + Gray ball
801109	TL01R	0.1ml-100ml	红色部件+灰色球型手握球 Red parts + Gray ball
801110	TL01G	0.1ml-100ml	绿色部件+灰色球型手握球 Green parts + Gray ball
801111	TL01C	0.1ml-100ml	灰色部件+灰色球型手握球 Gray parts + Gray ball
801112	TL02B	0.1ml-100ml	蓝色部件+黑色圆柱型手握球 Blue parts + Black cylinder
801113	TL02Y	0.1ml-100ml	黄色部件+黑色圆柱型手握球 Yellow parts + Black cylinder
801114	TL02R	0.1ml-100ml	红色部件+黑色圆柱型手握球 Red parts + Black cylinder
801115	TL02G	0.1ml-100ml	绿色部件+黑色圆柱型手握球 Green parts + Black cylinder
801116	TL02C	0.1ml-100ml	灰色部件+黑色圆柱型手握球 Gray parts + Black cylinder
801117	TL03B	0.1ml-100ml	蓝色部件+灰色圆柱型手握球 Blue parts + Gray cylinder
801118	TL03Y	0.1ml-100ml	黄色部件+灰色圆柱型手握球 Yellow parts + Gray cylinder
801119	TL03R	0.1ml-100ml	红色部件+灰色圆柱型手握球 Red parts + Gray cylinder
801120	TL03G	0.1ml-100ml	绿色部件+灰色圆柱型手握球 Green parts + Gray cylinder
801121	TL03C	0.1ml-100ml	灰色部件+灰色圆柱型手握球 Gray parts + Gray cylinder